

GTR-IA

Analog Tube Rotator


Quick Start Guide

What's in the box:

- Tube Rotator
- Power Adapter (w 3 each plug types)
- 1.5ml/2.0ml Tube Clips (12 each)
- 15ml Tube Clips (12 each)
- 50ml Tube Clips (12 each)
- Quick Start Guide

INITIAL SETUP

The tube holder plate is permanently mounted to the base unit. The tube holder plate has multiple channels for mounting various size tube holder clips. The clips can be installed in any configuration depending on the desired tube formats.

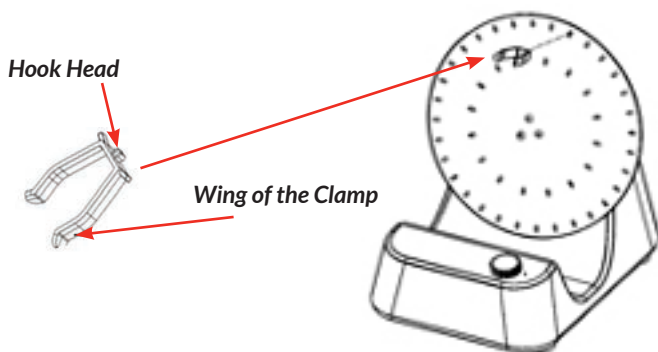


- Hold the two wings of the clamp, then place the hook head into the wide side of the desired channel and slide the hook head into the narrow end to lock in place. It is recommended that two tube clips be used for each larger volume (15mL and 50mL) tubes.

- Once the desired tube clips are installed, the GTR-HA will look similar to the image at right



Power Switch & Power Cord Socket Located on back of Unit

Speed Control Knob

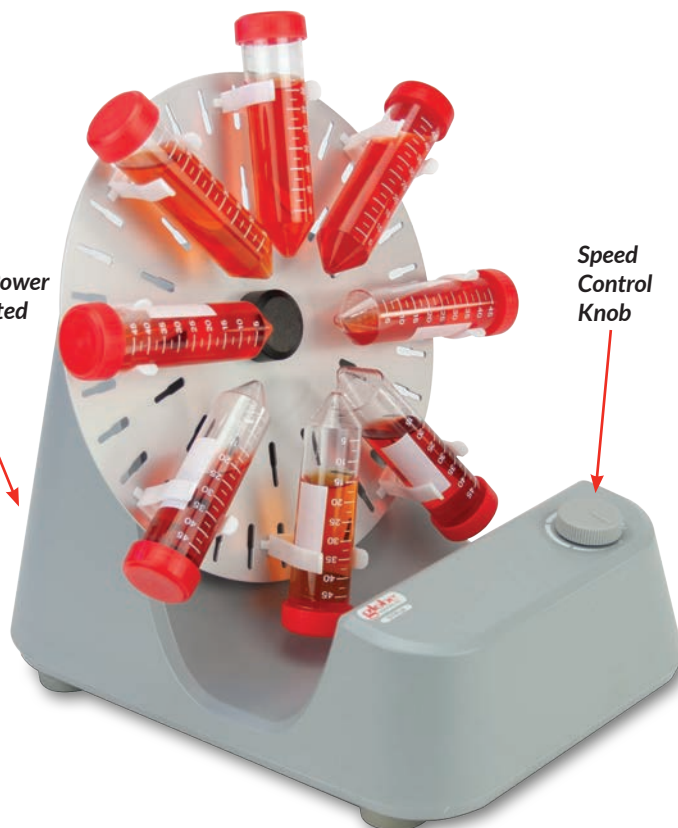

OPERATION

- Connect the power cable
- Install the tube holder clips as described at left
- Switch the equipment ON
- Rotate the speed control knob clockwise to set the desired speed and the actual rotation speed will adjust accordingly
- Rotate the speed control knob fully counterclockwise to turn off the rotation / mixing function

For the full User Manual, please visit www.globescientific.com

globe[®]
SCIENTIFIC INC.

Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400

Fax: 1 (201) 599-1406 • E-mail: mail@globescientific.com