


GTR-ID

Industrial Tube Rotator

Quick Start Guide


What's in the box:

- Quick Start Guide
- Tube Rotator
- 50mL Tube Holder Disk
- Power Cables (3 each)
- Set of (4 each) Screws, Flat Washers and Lock Washers
- Allen Wrench
- Extra Fuse

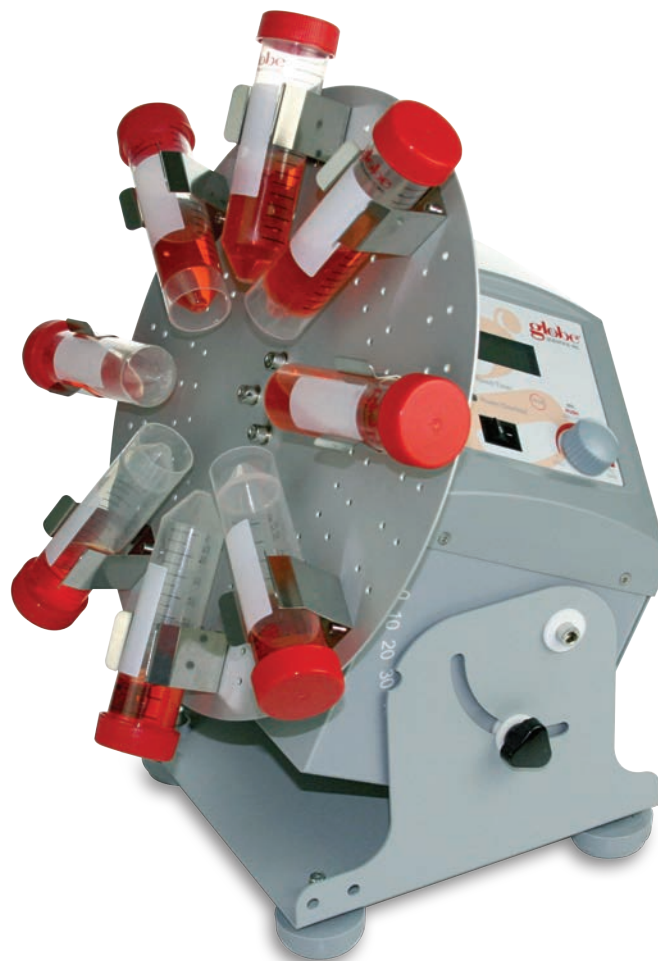

INITIAL SETUP

The 50ml tube holder disk comes standard with the GTR-ID, but it must be installed. The instructions below show how to install any available tube disk holder accessory:

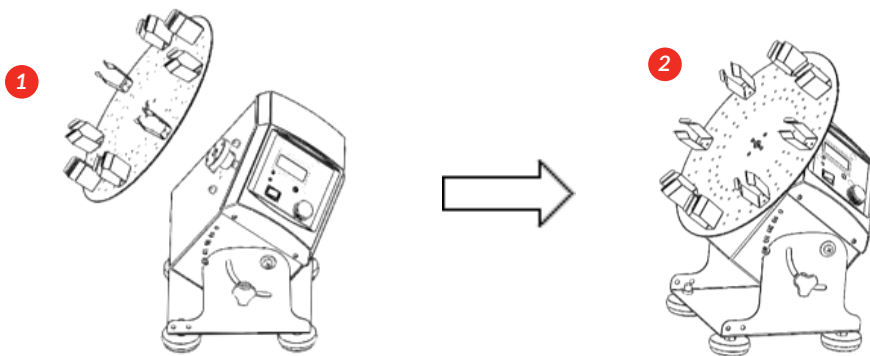

Secure the tube holder disk accessory with the four screws, lock washers and washers (installed in that order) using the included Allen wrench.

Continued ...

GTR-ID

Industrial Tube Rotator

Quick Start Guide

CONTROL PANEL / OPERATION



Speed / Timer Indicator
Green LED indicates the LCD screen currently displays speed. Red LED indicates the LCD screen currently displays time.

Power / Overload Indicator
Green LED indicates power on. Steady red LED indicates overload protection is enabled. Flashing red LED indicates overload and rotator operation is stopped.


On / Off Switch
Toggles power to the unit on and off.

LCD Display
Displays the working state, all set values, and any errors. (See table below.)

Mode Switch Button
Press to switch between speed and time.

Speed / Timer Control Knob
Rotate to set speed and time. Push to start/stop.

Display	Description
Set	Shown when setting a target value.
Err	Shown when an error occurs.
Display Area:	When Set is shown, the display shows the set value; When Set disappears, the display area shows the running value.



For the full User Manual, please visit www.globescientific.com



Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400

Fax: 1 (201) 599-1406 • E-mail: mail@globescientific.com